

Surface Evaporator

Made of opaque fused silica





Wissenschaftliche Apparaturen und Industrieanlagen AG Oerlikonerstrasse 88, CH-8057 Zürich Postfach 6962, CH-8050 Zürich Tel. 044 317 57 57 Fax 044 317 57 77 http://www.wisag.ch e-mail: info@wisag.ch

Surface Evaporator made of opaque fused silica

for rapid, economical evaporation and concentration of liquids and solutions

A helically electrical filament winding, covered by opaque fused/ transparent silica, with a big surface which is assembled in an opaque fused silica paraboloid dish with a grip, permits an indirect temperature rise heating of liquids by infra-red radiation.


In this way, rapid, economical evaporation or concentration of liquids is possible.

According to this principle, only the upper surface of the liquid to evaporate is heated very quickly as most of the radiation heat is absorbed already in low depth of penetration and, thus, it enables an economical, regular evaporation. Contrary to simmering evaporation, there is no substance loss because of splashing or super-boiling.

As surface evaporators are exclusively made of opaque fused/ transparent silicia, it is possible to make sensibly use of this material and its many advantages, as there are good IR-permeability, heat resistance, resistance to changes of temperature, extremely low electrical conductivity, exceptionally high chemical purity and resistance. These advantages of the material permit an economical evaporation of acids and other liquid chemical, except for hydrofluoric acid, phosphorus acid, and liquids with high alkali contents.

Electrical Safety

The opaque fused silicia encasing tube is a very good electrical insulator – protection against electric-shock hazard, optimal electrical safety by means of earthing of the inner tube. Surface evaporators meet the EG standard 89/336/EWG and EG standard 73/23/EWG.

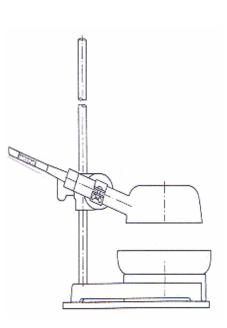

Surface	Order No.	Power	Dimensions				Net Weight	
Evaporator Type		Wattage	D	L	Н	d	approx. kg	
		Ū	mm	mm	mm	mm	1	1 The second
OV 140	9705011	700	140	120	55	27	1.0	i olt
OV 200	9705021	1000	200	190	70	40	1.4	
OV 300	9705031	1500	300	190	100	40	2.1	
Evaporation	Order	Volume	Dimensions				Net Weight	Н
basin	No.						• · · · · · · • · ·	+
ROTOSIL		ccm	D		Н		approx. kg	
flat version			mm		mm			
to OV 140	9705012	400	160		50		0.4	All surface evaporators are
to OV 200	9705022	1000	230		70		0.8	delivered with a silicone
to OV 300	9705032	4000	350		100		2.3	connection piece and a 1.5 m
Tripod	Order	Base	Dimensions		Net Weight	long cable with earthing		
	No.							
		Area	Н				approx. kg	contact-type plug of PVC.
		mm		mm				
to OV 140	9705017	330 x 250		520			3.5	They are designed for
								230 V ~.
to OV 200	9705017	330 x 250		520			3.5	
to OV 300	9705017	330 x 250		520			3.5	



Wissenschaftliche Apparaturen und Industrieanlagen AG Oerlikonerstrasse 88, CH-8057 Zürich Postfach 6962, CH-8050 Zürich Tel. 044 317 57 57 Fax 044 317 57 77 http://www.wisag.ch e-mail: info@wisag.ch